

15. ULTRASONIC CLEANER DEGAS SWEEP

Model	DR-DS13	DR-DS20	DR-DS30	DR-DS40	DR-DS45	DR-DS60	DR-DS100	DR-DS120	DR-DS150	DR-DS280
Volume Size (L)	1.3	2	3	4	4.5	6	10	12	15	28
Tank Internal Dimension (mm)	150 x 137 x 65	150 x 137 x 100	238 x 138 x 100	302 x 152 x 100	302 x 152 x 150	302 x 152 x 150	298 x 239 x 150	490 x 138 x 150	325 x 295 x 150	498 x 296 x 200
Operation Frequency	37kHz	37kHz	37kHz	37kHz	37kHz	37kHz	37kHz	37kHz	37kHz	37kHz
Ultrasonic Power (W)	80	80	160	160	160	240	320	400	480	640
Heating	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓
Degas	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓
Sweep	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓
Drainage Valve	NO	NO	O	O	O	O	✓	✓	✓	✓
Carrying Handle	NO	NO	✓	✓	✓	✓	✓	✓	✓	✓
Lid	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓
Unit Weight (Kg)	2.2	2.3	3.2	5	4.5	5	7	7	13	19
Unit / Carton	4	4	4	2	2	2	1	1	1	1
Carton Size (cm)	53 x 53 x 35	53 x 53 x 35	49 x 34 x 60	43 x 57 x 34	49 x 34 x 39	57 x 42 x 39	47 x 37 x 43	65 x 25 x 47	47 x 42 x 47	65 x 42 x 49
Carton Weight (Kg)	12	13	17	11	10	13	9	9	15	21

✓ : Included ○ : Optional NO : Not Available

Features Illustration:

DR-DS30



7) Window shows the minute just set or minute remained in ultrasonic cleaning process

6) Actual °C/°F window shows the current temperature of cleaning fluid inside the tank

8) Timer setting

5) Heating setting

4) Sweep button

3) Degas button

2) START/PAUSE & Memory Setting

1) Power ON/OFF

16. OPERATION INSTRUCTIONS:

- 1) Check the specification table for the correct operating requirements (located on the rear of the unit).
- 2) Before connecting power, put the unit on a stable and flat working surface. Add water to tank, suitable to completely immerse the items to be cleaned. Using stainless steel basket is highly recommended.

! Never operate the unit without water in the cleaning tank. Always fill the unit above 2/3 marker line indicated on the inside of the tank. Failing to do so, DAMAGE MAY OCCUR.