

FD Series Vertical Autoclave


Product Features

1. Simple and easily understandable control panel to show the sterilization temperature, sterilization time, drying time, press separate keys to set the temperature and time makes the operation easier
2. Each sterilization stage is displayed with LED lights, making the sterilization process clear at a glance.
3. Microcomputer intelligent control system ,achieves fully automatic control of the sterilization process.
4. Pulse method to completely purge out cold air.(**E/R only**)
5. System automatically monitors the discharge of cold air to ensure a pure steam sterilization environment and the best sterilization effect.
6. Up to 20 error information records can be stored, facilitating instrument management.
7. Built in water tank for internal discharging of steam, effectively absorbing water vapor from exhaust process.(**E/R only**) Fully automatic drying function or semi-automatic drying function choice(**R only**)
8. Fully automatic drying: Equipped with an air pump and a 0.2um air filter,it automatically starts drying after sterilization. After sterilization, clean air will be put in to replace the wet air inside, so as to improve drying efficiency.(**Fully automatic drying is an option for R only**)



Safety Features

- **Lid closing checking:** Machine can not start if lid is not locking well,the self-induction interlock device can avoid risk of misuse.
- **Unique anti-scald cover:** Nice look and avoid causing scald.
- **Triple protection for lack of water:** The bottom of the sterilization chamber is equipped with three different dry burning protection devices: liquid expansion type, copper temperature sensing type, and electric ion type (water level sensor), to avoid misjudgment caused by a single method.
- **Double protection for over pressure:** Mechanical pressure gauge and electric sensing system, once pressure is abnormal, system will cut power, release the pressure and send out alarm.
- **Auto trouble shooting:** System will monitor the working status, and will cut the power, send out alarm, and report the error information once any abnormal situation detected.
- **Safety test function:** Regular testing of system security and safety protection measures is possible.
- **Real time monitor:** System will monitor the lid locking, pressure and temperature status on a real time basis to guarantee the safe and successful work of the autoclave.
- **Electrical protection:** Equipped with a highly sensitive electrical protection system, effectively providing leakage protection, overcurrent and short-circuit protection.

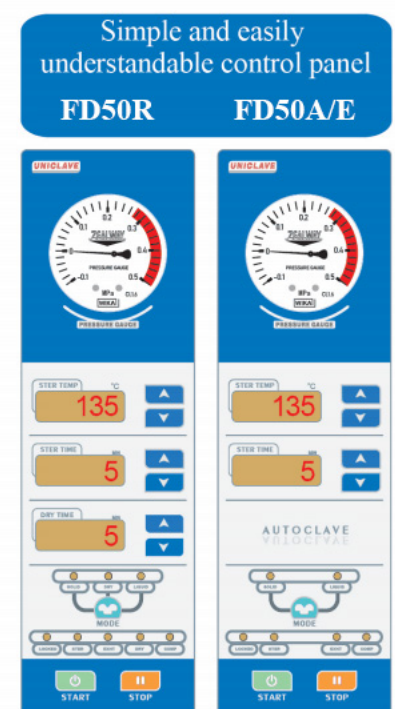

Specification

| Model | Basic | FD29A | FD50A | FD80A | FD100A |
|------------------------------|---|---|---------------|------------------------|---------------|
| | Pulse | FD29E | FD50E | FD80E | FD100E |
| | Drying | | FD50R | FD80R | FD100R |
| Difference | FD-A: Fully automatic FD-E: FD-A+pulse+inner steam tank+waste water tank FD-S: FD-E+semi automatic drying function, if need automatic drying need to add 530USD | | | | |
| Capacity(L) | | 29 | 50 | 80 | 100 |
| Dimension(L*W*H)(mm) | | 580x490x960 | 580x490x 1085 | 730x625x1100 | 730x625x1200 |
| Chamber dimension(Dia*H)(mm) | | Φ 320x378 | Φ 320 x725 | Φ 400 x700 | Φ 400 x 800 |
| Power requirements | | 220V±10% 16A 50HZ/60HZ | | 220Y±10% 32A 50HZ/60HZ | |
| Rated power(W) | | 2600W | 2600W | 4600W | 4600W |
| Sus304 basket size(pcs) | | (Φ 298x260)x1 | (Φ 298x195)x3 | (Φ 360x275)x2 | (Φ 360x215)x3 |
| Chamber material | | SUS304 stainless steel | | | |
| Sterilization temperature | | 105°C~135°C | | | |
| Sterilization time range | | 1~300min | | | |
| Pressure gauge range | | -0.1MPa~0.5MPa | | | |
| Pulse times | | 0~9 times(E/R only) | | | |
| Exhaust temperature range | | 73°C~104°C(E/R only,liquid mode)105°C~130°C(Solid mode) | | | |
| Safety valve set pressure | | 0.29MPa | | | |
| Drying | | Drying temperature:80°C~160°C Drying time:1~300 min | | | |
| Sterilization mode | | Solid mode Liquid mode Sterilization+ drying(R only) | | | |
| Controller | | Microcomputer control system | | | |
| Safety device | | Lid closing checking, over temperature protection system,dry scorch protection , over pressure protection, safety valve, over current/short circuit/electric leakage protection,scald protection, automatic failure-inspection system | | | |

Note:FD-A has liquid mode, but can not preset the exhaust temperature for liquid mode.

Spare parts option



Micro printer
To print temp.,time
and F0 value



Digital pressure sensor



Tray



Stainless steel bucket



IQ/OQ/PQ
(3Q)adapter joint



Adjust foot

Printing set:printer+digilal
pressure sensor:can print temp. and
pressure in data and flow chart way